

RheBuild® Kits

RheBuild Kits are available for all Rheodyne® brand products. Included in each individualized RheBuild Kit are all parts, tools, and instructions to maintain precision performance of your particular product. RheBuild Kits eliminate individual part ordering.



APPLICATION NOTE

How to Avoid Pressure Transients

Air in the sample loop can cause an instantaneous system pressure drop that eventually returns to a normal level. Air causes the pressure to drop when the injector moves from the LOAD to the INJECT position. When large sample loops ($\geq 100 \mu\text{L}$) are partially loaded, air present in the needle port tube is pushed into the sample loop (see Figure 1). Air can also enter the sample loop from siphoning which occurs when the vent line is higher than the injection port. In either case, upon injection, the system pressure collapses the air bubble, causing pressure to drop momentarily.

A pressure drop in the system caused by air results in changes in retention time, artifact peaks, and affects column performance.

Avoid pressure drops by removing the air in the needle port tube. Do this by flushing about 1 mL of mobile phase with a luer syringe with needle port cleaner. Keep the needle port tube filled with mobile phase by occasional flushing. Adjust the vent line(s) so the outlet is at the same horizontal level as the needle port (see Figure 2). For additional injection troubleshooting, refer to the Rheodyne Troubleshooting Guide for HPLC Injection Problems. You may download the Guide from the IDEX Health & Science web site: www.idex-hs.com under Support.

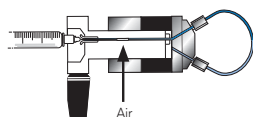


Figure 1 Air present in the needle port tube is pushed by the syringe during loading into the sample loop

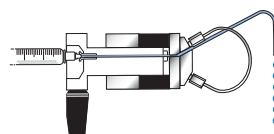


Figure 2 Pathway of the flushing mobile phase using the Needle Port Cleaner, Part # 7125-054 (see page 143) when the injector is in INJECT

Part No. Description

RHEBUILD KITS FOR MX SERIES II™ VALVES

7150-999	RheBuild Kit for MXT715-000 (includes 2 rotor seals)
7152-999	RheBuild Kit for MXT715-102 (includes 2 rotor seals)
7155-999	RheBuild Kit for MXT715-105 (includes 2 rotor seals)
7920-999	RheBuild Kit for MXP7920-000 and MXP7900-000
7960-999	RheBuild Kit for MXP9960-000 (includes rotor seal and stator face seal)
7961-999	RheBuild Kit for MXP7960-000
7970-999	RheBuild Kit for MXP7970-000
79801-999	RheBuild Kit for MXP7980-000
79861-999	RheBuild Kit for MXP7986-000
7900-999	RheBuild Kit for MXP9900-000 (includes rotor seal and stator face seal)

RHEBUILD KITS FOR MANUAL VALVES

3725-999	RheBuild Kit for models 3725, 3725i, 3725-038, 3735i-038
7010-996	Conversion Kit including Stator Face Assembly for model 7010
7010-997	RheBuild Kit including Stator for model 7010
7010-998	RheBuild Kit, pH Upgrade Kit for model 7000
7010-999	RheBuild Kit for model 7010 and 7010-type Valves
7125-999	RheBuild Kit for models 7125 and 7126
7125Ti-999	RheBuild Kit for model 7125-081
7410-999	RheBuild Kit for model 7410
7520-999	RheBuild Kit for models 7520 and 7526 (includes inlet stator and seal)
7725-999	RheBuild Kit for models 7725 and 7725i
7788-999	RheBuild Kit for model 7725i-188
8125-999	RheBuild Kit for models 8125 and 8126
9010-999	RheBuild Kit for model 9010
9125-999	RheBuild Kit for models 9125 and 9126
9725-999	RheBuild Kit for models 9725 and 9725i; 7725(i) pH upgrade kit

RHEBUILD KITS FOR MX SERIES I™ VALVES

7900-999	RheBuild Kit for models MX7900-000, MX7925-000, MX9900-000, MX9925-000
7960-999	RheBuild Kit for model MX7960-000
7980-999	RheBuild Kit for model MX7980-000
7984-999	RheBuild Kit for model MX7984-000
7986-999	RheBuild Kit for model MX7986-000

RHEBUILD KITS FOR LABPRO™ & EV AUTOMATED FLUIDIC INSTRUMENTS

1006-999	RheBuild Kit for model PR/EV100-106
5001-999	RheBuild Kit for models PR/EV500-101 and PR/EV550-101
5100-999	RheBuild Kit for models PR/EV500-100 and PR/EV550-100
5104-999	RheBuild Kit for models PR/EV500-104 and PR/EV550-104
7004-999	RheBuild Kit for models PR/EV700-104 and PR/EV750-104
7112-999	RheBuild Kit for models PR/EV700-112 and PR/EV750-112
7501-999	RheBuild Kit for models PR/EV700-100 and PR/EV750-100
7502-999	RheBuild Kit for models PR/EV700-102 and PR/EV750-102
7507-999	RheBuild Kit for models PR/EV700-107 and PR/EV750-107
7531-999	RheBuild Kit for models PR703-100 and PR753-100